


HOSOKAWA MICRON GLOVE TESTER

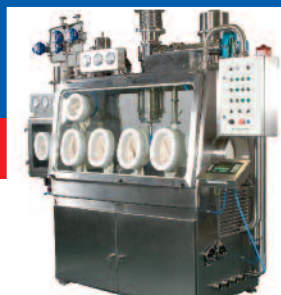

Hosokawa Micron's new automated glove integrity tester provides a quick and easy way to test isolator fitted pharmaceutical gloves and gauntlets.

This fully automated, easy to use glove tester is essential to check for damage to gloves and gauntlets, not apparent to the naked eye, which may put operators handling potentially harmful materials at risk.

- Clear and simple glove integrity verification
- Even detects holes less than 100µm
- Simple to use. Two button operation
- Fully automated measurement
- Wireless, lightweight, hand held device
- Automatic glove filling
- No additional air supply required
- Ready for use straight from the case
- Complete with battery, battery charging stand and case



With preconfigured parameters tailored to your glove type and detection requirements the Hosokawa Micron Glove Tester is quick and simple to use either before each procedure or as part of scheduled integrity tests.

Operation:

Clip battery into place on hand set. Place handset into glove port. Press START.
Handset seal inflates to fit gloveport, clamping handset in place.
Air is used to inflate gauntlet.
A digital display confirms pressure and pass/fail.



HOSOKAWA MICRON LTD

Rivington Road, Runcorn WA7 3DS Tel: +44 (0)1928 755100 • e.mail: info@hmluk.hosokawa.com • www: hosokawa.co.uk

PROCESS TECHNOLOGIES FOR TOMORROWSM